

VWR-214B2 Analytical and Precision Balance

Note: closing the doors will mitigate inaccuracies resulting from air draft and instrument imbalance

- Press the “ON” button to turn the instrument on
- Slide open the door and gently place an empty weigh tray onto the weigh plate
- Close the door and tare the balance by pressing the “→T←” button on the terminal
- Open the door and carefully load the sample onto the weigh tray using a spatula
- Close the door and receive the final weight
- Clean any excess sample using a wipe
- Turn off by pressing the “OFF” button

