

Mettler Toledo WXTS3DU Microbalance

Important notes:

- The microbalance is highly sensitive so do not press on the weigh plate or overload the weigh plate
- For greatest accuracy do not exceed 10g
- Closing the doors will mitigate inaccuracies resulting from air draft and instrument imbalance
- Avoid suddenly impacting the weigh plate

Operation:

1. Press the power button on the terminal to turn the instrument on and then press the “→0←” button
2. Slide open the door and gently place an empty weigh tray onto the weigh plate
3. Close the door and tare the balance by pressing the “→T←” button on the terminal
4. Open the door and carefully load the sample onto the weigh tray using a spatula
5. Close the door and receive the final weight
6. Clean any excess sample using a wipe
7. Turn off by pressing the power button

